User Manual

© 2004, Xcitex, Inc., All rights reserved.


Overview

The Professional Analyst, Flight Path X-Ray Edition software can provide accurate 3-D measurements of points and angles using simultaneously acquired images from a pair of X-Ray systems. This document describes the application interface, the calibration and measurement procedure, and the data export capabilities of the software.

For more information on the technical background of this software and a discussion of the accuracy and sources of error, please refer to the *Professional Analyst, Flight Path Edition, Brief Technical Overview* document.

Application


The application has three main document types: Calibration, Measurement, and Event Summary. The Calibration document uses images of a calibration fixture taken by a pair of X-Ray systems and computes a calibration for each pair. The resulting calibration pair is used by the Measurement document to do precise measurements on a pair of images from an event. A sequence of Measurement documents is used to construct an Event Summary document that compares the results from each of the Measurement documents to produce a summary of the event.

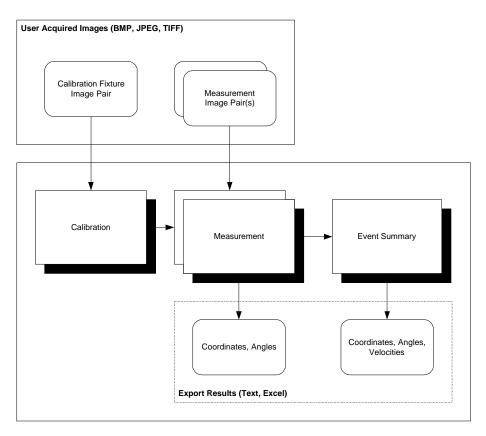

Each of these document types will be explained in more detail in the later sections.

Application Windows

The main application window has three primary sections: the main desktop area where documents appear, the Graph windows, typically docked on the right, and the Points List, typically docked on the bottom.

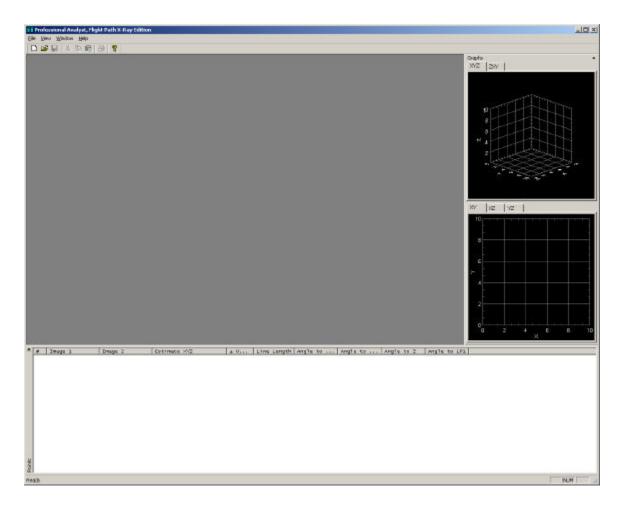

The Graphs dialog contains a set of 3D graphs and a set of 2D graphs. The 3D graph can be rotated by clicking and dragging the left mouse button. To pan the 3D and 2D graphs, hold the **Shift** key while clicking and dragging the left mouse button. To zoom the 3D and 2D graphs, hold the **Alt** key while clicking and dragging the left mouse button. If your mouse is equipped with a wheel button, the wheel can also be used to zoom the graph windows. Additional options are available in the context menu by clicking the right mouse button over the graphs.

The Points List provides information about all selected points for the Calibration and Measurement Documents and it provides information about the Measurement Line Pairs for the Event Summary Documents.

Document windows typically have controls on the right and images on the left. The images can be panned and zoomed using the same controls as the graphs. **Shift** + left mouse button for panning, and **Alt** + left mouse button for zooming. The mouse wheel also controls zooming. When using the mouse wheel to zoom, the image will zoom in/out on the currently selected point. This allows the user to very quickly zoom in and out on a particular feature. Click where you would like to zoom, then use the mouse wheel. The application will automatically center the current selected point.

Document windows will have a toolbar across the top of each window. The document toolbar typically contains some or all of the following controls. Many of the items below have separate settings for the top and bottom views. Click on the view that you wish to change the settings for and then click on the desired toolbar button. A green box in the upper right corner of each view indicates which view is currently active.

Toggle Side Panel

Switch the side panel between a small toolbar version and the full size version.

📁 Image Settings

Adjust image settings such as brightness, contrast, gamma, etc.. This control has separate settings for the top and bottom views.

Locate Point Settings

Adjust parameters for the automatic location of points for Calibration Documents. This feature will be discussed later in this document. This control has separate settings for the top and bottom views.

Define World Reference Frame

Define the transformation to the world reference frame for Calibration Documents. This feature will be discussed later in this document.

Enable Midpoint Finder

Enable the midpoint finder tool. This tool allows you to click and drag two endpoints of a line in the image. The program will draw a marker that indicates the center of the line. The program will also draw a circle through the two endpoints and centered on the center of the line. To hide the midpoint finder tool, click anywhere in the image, but do not drag.

Enable Reticle

Show or hide the reticle (cross-hairs) in the view. This control has separate settings for the top and bottom views.

Enable Rulers

Show or hide rulers along the edges of the view. This control has separate settings for the top and bottom views.

Show Point Labels

Show point labels (numbers) next to each point in the view. This control has separate settings for the top and bottom views.

G Flip Vertical

Flip the image vertically. This control has separate settings for the top and bottom views.

Flip Horizontal

Flip the image horizontally. This control has separate settings for the top and bottom views.

11 View Actual Size

Resize the displayed image so that it is shown at the actual resolution of the image. This control has separate settings for the top and bottom view.

🗵 Fit in Window

Resize the displayed image so that the complete image is shown in the view. This control has separate settings for the top and bottom view.

+ Pan Mode

Set the mouse mode to pan mode so that all left click and drag mouse operations will pan the image. This control has separate settings for the top and bottom views.

Zoom Mode

Set the mouse mode to zoom mode so that all left click and drag mouse operations will zoom the image. This control has separate settings for the top and bottom views.

Typical Test Scenario

A typical test scenario has multiple steps. The first step is calibration. During calibration, you prepare the X-Ray systems to be able to perform precise measurements by determining properties of the X-Ray systems. This is typically done by capturing images of a calibration fixture and then analyzing these images to produce a calibration file. The next step is to run the actual projectile tests to conduct measurements. During measurement, images of the projectile are captured by the calibrated X-Ray systems. These measurement images, combined with the calibration file, are then used to produce 3D measurements of points on the projectile.

The calibration step only needs to be performed as necessary. If none of the X-Ray systems have been moved or changed since the last calibration, then the previous calibration should be satisfactory. The uncertainty in the measurements can be used to determine if a new calibration is required. If the uncertainty exceeds a desired limit, then a new calibration must be performed.


If there are multiple X-Ray stations located along the firing path, then calibration and measurement must be performed for each station. The measurements from each of the X-Ray stations are then combined to produce an event summary. An event summary is a collection of measurements from multiple camera or X-Ray stations for a single firing event. The event summary computes additional information by comparing the results from each of the measurements, such as velocity and angular velocity.

Calibration

In order to perform accurate measurements using the X-Ray system, a calibration must first be performed. This calibration computes necessary information about the properties

and location of each X-Ray system. If any of the systems are moved or changed after calibration, then a new calibration must be performed.


Calibration is performed by capturing images of special *calibration fixtures* that have fiducial markers at precisely known locations. These images can then be analyzed and the properties of the X-Ray systems can be determined.

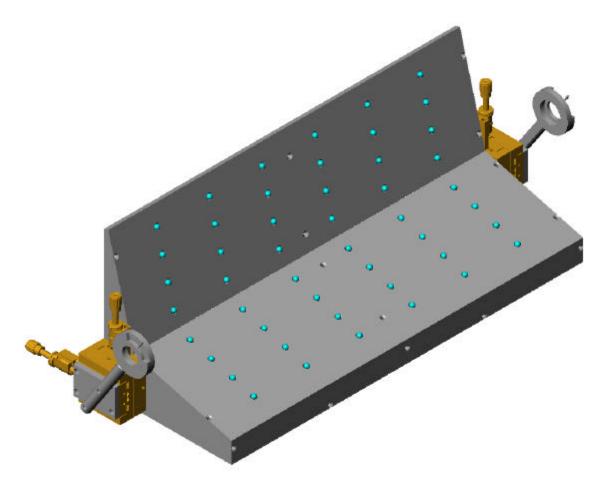

A typical calibration procedure for an event with multiple stations:

- 1. A laser should be used to ensure alignment of each calibration fixture relative to the firing path. Place the laser next to the gun barrel and shine it down the expected path of the projectile.
- 2. Place the calibration fixture in the center of the field of view of each X-Ray system, roughly aligned to the path of the projectile (within ±6-7mm). The entire calibration fixture (including irises) should be completely visible in both X-Ray views.
- 3. Adjust the position of the irises on the ends of the calibration fixture using the fine adjustment screws so that the laser passes through the exact center of both irises when they are closed. If the laser cannot be centered on the irises, then shift the whole calibration fixture so that the laser can be centered in the irises.

- 4. Measure the distance from the gun barrel to the first iris on each calibration fixture. This will be the *Z Coordinate of the First Point* for each fixture. This value is used to help define the world reference frame.
- 5. Measure the angle of each calibration fixture in the direction perpendicular to the path of the projectile. This will be the *Base Angle* for each fixture. This value is used to help define the world reference frame.
- 6. Capture images of each calibration fixture from each X-Ray system.

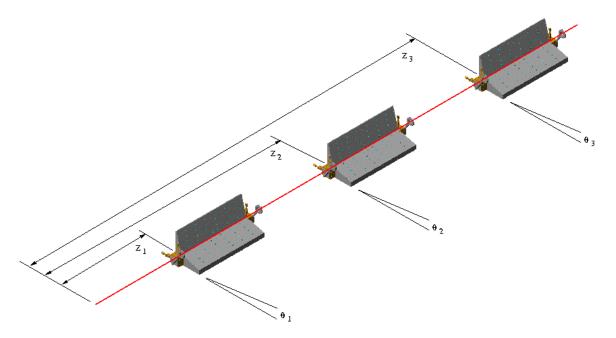


The resulting images are then used to create Calibration Documents for each X-Ray pair.

A Calibration Document is created by selecting **File** \rightarrow **New**.



Then select **Calibration**. An empty calibration document will appear.

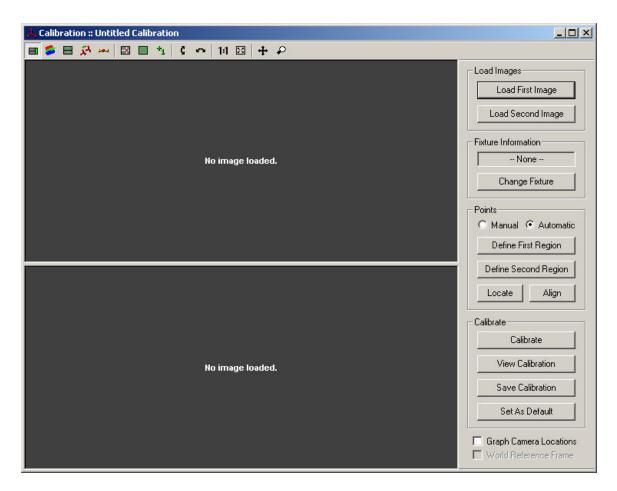

All controls are located on the top toolbar and on the left panel of the Calibration Document window. The toolbar controls were described earlier in this document.

The standard procedure for calibrating X-Ray systems is as follows:

- 1. Load the first calibration image by clicking on the **Load First Image** button and select the correct image file (BMP, TIFF, or JPEG). This image should be the top view looking down upon the calibration fixture.
- 2. Load the second calibration image by clicking on the **Load Second Image** button and select the correct image file. This image should be the side view looking horizontally at the calibration fixture.
- 3. Load the fixture information file for the specific calibration fixture that is being used. The fixture file (*.fixt) should have been supplied with each calibration fixture. The fixture file contains information about the location of the fixture points and the default units of measure.
- 4. Now the fixture points need to be identified in the top and bottom images. There are multiple options for locating the fixture points. It can be done manually or automatically.

To locate points automatically:

- Select **Automatic** in the Points section of the right panel.
- Define a region around all the fixture points in the top image by clicking on **Define First Region**. Then click and drag the mouse to draw a rectangle around all the fixture points in the top image. If you make a mistake, click on **Define First Region** again and redraw the rectangle.
- Define a region around all the fixture points in the bottom image by clicking on **Define Second Region**. Then click and drag the mouse to draw a rectangle around all the fixture points in the bottom image.
- L Calibration :: xraytest1and5.calb - 0 × 🔳 🎓 🖶 📯 🚧 🔝 🔳 🐴 🕻 🗛 111 🖾 🕂 🖓 RETICLE = [1564.0, 948.0] Load Images Load First Image + +++++ Load Second Image Fixture Information xraytest1and5.calb + Change Fixture ÷ Points C Manual 💿 Automatic + ÷ Define First Region Define Second Region RETICLE = [118 8.0, 1241.0 ÷ Г + Locate Align Calibrate + + Calibrate ÷. + View Calibration Save Calibration ÷ + Set As Default + 🔲 Graph Camera Locations 🗖 World Reference Frame
- Click on **Locate** to automatically locate fixture points in both images.

- If the automatic location of points fails, there are parameters that can be adjusted to change the behavior of the automatic location scheme.
 - i. Click on the **Locate Point Settings** icon on the toolbar, or rightclick on the top or bottom image and select **Locate Point Settings** from the drop down menu.

C Black Points 🛛 🧿	White Points
Intensity Threshold	64
Minimum Pixel Count	100
Maximum Pixel Count	5000
Maximum Eccentricity	5
 Use Image Settings Use Edge Detection 	

There are different settings for the top and bottom images, so if necessary, change both settings separately.

 ii. In addition, improving the contrast of the fixture points can also help in automatically locating them. You can change the image settings for the top and bottom images by clicking on the Image Settings icon on the toolbar, or right-click on the image and select Image Settings from the drop down menu.

B&W Image Settings	
Driekhnana /	Reset
Brightness	
Contrast 🚽	Load From Fil
Gamma 🖣	1.0 Court To Th
Exp./Log. 🗸 📕	67 Save To File
Reverse	and a second

iii. If the image that has been loaded is a 16-bit TIFF file, you can also select which bits of the TIFF file to display. Right-click on each image and select **TIFF Bit Selection**.

Image Settings Locate Point Settings Define World Reference Frame Epoble Midpaint Finder	
Enable Midpoint Finder Enable Reticle Enable Rulers	_
Show Labels Flip Horizontal Flip Vertical	_
View Actual Fit in Window	_
Pan Mode Zoom Mode	
TIFF Bit Selection	Show Bits 1-8 Show Bits 2-9 Show Bits 3-10 Show Bits 4-11 Show Bits 5-12 Show Bits 6-13 Show Bits 7-14 Show Bits 8-15 Show Bits 9-16
	✓ With Saturation

To locate points manually:

- Select **Manual** in the Points section of the right panel.
- Click on a point in the top image corresponding to a fixture point.
- Click on a point in the bottom image corresponding to the same fixture point.
- Click on the **Add Point** button.
- If you wish to adjust the coordinates of a point, select the point in the Points list. Click on the correct points in the top and bottom images. Then click on the **Update Point** button.

	ensional Analysis, Flig na Wildon Bilo	ht Path X-Ray Edition						LO X
		91 8 1						
	Collection	r martest Landiticals	-					0-apts - XVZ Z/V
				P		10.995	ALC: U	
	REDCLE - 1956	44,9184				-Load	(mage)	
						- 1	Lost fittings	and the state
						- 1 <u>1</u> 1	and Second Image	s
							z biloandian	
					•		Harlen ficab	
					•		Change Rinue	
						- Park	l Hanad (? Automatic	
					•	100	Define Fact Region	
						- In	dire Second Region	W NZ NZ
	METHOLE - 1918	•			•	7 4	ecele Align	
						- Callor		
					•		Calibrate	11
					•		View Talibusion	324
					•		Save Collection	1
					•		SechaDalad	Fui
						E B	raph Cemera Locations	12
		-	÷	-		E.5	ord Patriction Frank	
								4 28 16 04 18 2 X[echec]
	Imoge 1	Drage 2	Calibration 202	Estima	2012	Error		
1	1315.5, 555.1	490.2, 822.7 939.9, 824.3	2.23, 4.45, 3.25, 4.65,	0.00 1.28, 3.00 1.25.	4.60, D.DF	0.01		
2	1035.0, 543.4	1367.2, 625.6 1845.1, 927.1	3.25, 4.65, 3.25, 4.65,	5:00 1:24, 9.00 1.23,	4.64, 6.00	0.01		
5	824.8, d18.4 1324.8, 622.6	488.3, 1001.8 938.5, 1004.0	8.80, 8.44, 8.90, 3.44,	0.03 0.82, 3.00 0.92,	E.45, 0.02	0.01		
2	1928.7, 623.9	1399, 1006	8.91, 3.44,	6.00 0.82,	2.45, 5.99	0.01		
2	3331.0, 618.0 841.6, 075.5	18+7, 1008 480.1, 1185.2	8.80, 2.23,	P.00 0.82, 0.00 0.61,	1.25, -0.00	0.01		
10	1222.5, 676.7	907.0, 1107.4 1391, 1191	8,60, 2.17,	3,00 0,61, e.00 0.ec,		0.01		
12	2328-1. 685.0	1841, 1150	8.80, 2.23,	9,00 0,40.	2.22. 8.99	0.01		
14	057.9, 772.9 1228.9, 727.1	47710, 1209.2 917.9, 1274.5	8.27, 1.07, 1.27, 1.02,	0.00 0.28, 1.00 0.28,	1.01, 2.99	0.02		
14	1802.2, 740.1 2315.1, 740.2	1540, 1301	8.27, 1.08, 8.27, 2.03,	z.03 0.27, 9,00 0,29,		0.01		
17 17 17 18	069.7, 948.5 1246.2, 948.5	458.9, 1572.0	-1.01, -0.17, -1.01, -0.17,	0.00 -1.15.	-3.26, -0.07	0.02 0.01		
alis	1425.9. 951.T		-1.mn.37.	6.01 -1.14		0.07		1
Reath								FA.M //

- To remove a point, select the points to remove in the Points list. Then click on the **Remove** button.
- After you have set all the fixture points, if you wish for them to be automatically aligned to a standard grid pattern used by typical calibration fixtures, then click on the **Align** button.
- 5. Once all fixture points have been located, click on the **Calibrate** button.

Professi	onal Analyst, Flig	ht Path	X-Ray Edition	x
(j)	Calibration Comp	leted:		
4	Mean Error	=	0.0254 in	
	Std. Deviation	=	0.0023 in	
	Min Error	i ₩	0.0106 in	
	Max Error	÷	0.0472 in	


After a pair of X-Ray systems have been calibrated, you can view the results of the calibration by selecting **View Calibration**.

irst X-Ray System		Second X-Ray System
Location		Location
-0.8630 60.204	3 5.9397	-44.4255 1.9811 5.1861
Rotation		Rotation
0.0044 0.007	9 1.0000	-0.0107 0.0049 0.9999
-0.9623 -0.271	8 0.0064	-0.2885 -0.9575 0.0016
0.2718 -0.962	3 0.0064	0.9574 -0.2885 0.0116
Translation		Translation
-6.4115 15.493	2 58.1329	-5.6690 -10.9283 43.0451
Center of Projection		Center of Projection
1893.05 -1597.7	'9	1339.85 3114.66
Axis Scale Factors		Axis Scale Factors
9258.75 8930.3	34	6489.71 6257.02
alibration Statistics		
Mean Error	0.0254	Standard Deviation 0.0023
Minimum Error	0.0106	Maximum Error 0.0472
		nches OK

After the X-Ray systems have been calibrated, for every point that you click on in the top image, an epipolar line will be drawn in the bottom image and vice versa. The epipolar line indicates which points in the bottom image could correspond with the point that you have clicked on in the top image.

The current point in the top image is shown with a red plus symbol. The corresponding epipolar line in the bottom image is shown as a red line. The current point in the bottom image is shown with a blue plus symbol. The corresponding epipolar line in the top image is shown with a blue line.

The red plus and the blue line and the blue plus and the red line may not exactly intersect. The red and blue lines represent images of two rays coming from each X-Ray system. Since they do not exactly cross, the midpoint of the shortest distance between the two lines is shown with a yellow plus symbol in both the top and bottom views. The distance from the yellow plus to the red and blue lines is the uncertainty in the position estimate. Information along the bottom of each view indicates the 3D coordinates of the midpoint and the uncertainty associated with the 3D coordinates.

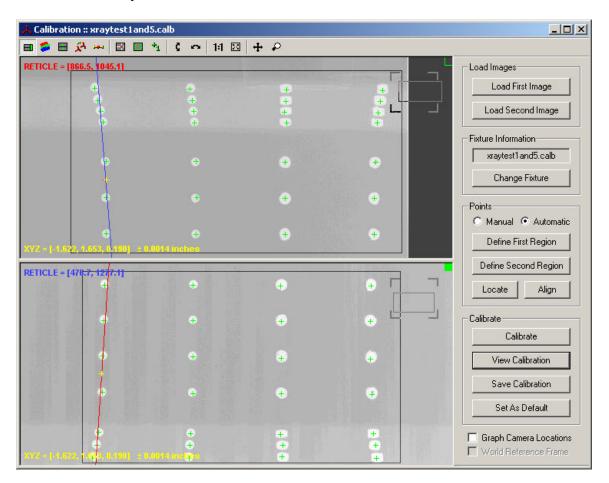

When you are satisfied with the calibration results, you can save the calibration to a file using the **Save Calibration** button. This will save all the calibration information to a calibration file (*.calb) that can be used by the Measurement document to perform measurements on other images taken by the same X-Ray systems.

If you would like the same calibration to be used for all newly created Measurement documents, then select the **Set As Default** button after the calibration has been saved to a file.

World Reference Frame

The coordinate frame of the calibration is the coordinate frame that is used by the calibration fixture file that was loaded. All the coordinates in the fixture file are referenced from some point of origin on the fixture. In order to compare information from one pair of X-Ray systems to another pair of cameras or X-Ray systems, the coordinates need to be referenced to some mutual world reference frame. All coordinates

can then be transformed from the local calibration fixture reference frame to the world reference frame.

A tool is provided to help define the transformation between the calibration fixture and world reference frame. You can access this tool by selecting the **Define World Reference Frame** button on the toolbar, or by right-clicking in the image and selecting **Define World Reference Frame**.

	anslation (2	<w *="" +<="" =="" r="" th="" xc=""><th>IW]</th></w>	IW]
Rotation (R)			
	-0.069754	Contraction of the second	
0.0697946	0.997546	-0.005475	
-0.006773	0.0059630	0.999959	
Translation (Tw	d.		Units
0	0	50	Inches 🔻
-1.69946	2.39674	-8.09993	Get Current
1		1	
Second Point			
Second Point -1.91497	2.56127	21.9444	Get Current
	line stateme	21.9444 Units	Get Current
-1.91497	line stateme		Get Current
-1.91497 Z Coordinate ol	f First Point	Units Inches	Get Current

There are two methods for entering information about the world reference frame. You can directly type in the rotation and translation matrix used to transform calibration coordinates to world coordinates. Or you can select two points that lie on the world reference frame Z-axis and supply the Z coordinate of the first point and the base angle of the fixture.

For the first method, if the rotation matrix and translation vector are known, simply enter them into the appropriate boxes and hit **OK**.

For the second method, the standard calibration fixture that was designed to be used with this software should be equipped with two adjustable irises, one on each side of the fixture. The calibration fixture should have been positioned so that the irises are aligned with the flight path and are visible in both the top and bottom views. To define the world reference frame:

- 1. The system should have been calibrated already, as described earlier.
- 2. Click on the center of the left iris in the top and bottom views of the calibration document window. Check the uncertainty in the coordinates to make sure that a suitable point has been chosen.
- 3. Click on **Get Current** for the **First Point** in the Define World Reference Frame dialog. An orange box should appear in the calibration document window to indicate the first point.
- 4. Click on the center of the right iris in the top and bottom views of the calibration document window. Check the uncertainty in the coordinates to make sure that a suitable point has been chosen.
- 5. Click on **Get Current** for the **Second Point** in the Define World Reference Frame dialog. An orange box should appear in the calibration document window to indicate the second point and a dotted line should be drawn between the first and second points.
- 6. Enter in the **Z** Coordinate of the First Point and the units of the Z coordinate. This should have been measured when the calibration fixtures were being aligned.
- 7. Enter in the **Base Angle** in degrees. This should have been measured when the calibration fixtures were being aligned.
- 8. Hit **Apply** and the equivalent rotation and translation will be computed and applied.


Once the world reference frame has been defined, the coordinates of all points can be displayed in either the local calibration reference frame or the world reference frame. You can select which frame to display in by checking the **World Reference Frame** checkbox in the side control panel. When displaying coordinates in the world reference frame, the graphs will label the axes as Xw, Yw, and Zw.

The world reference frame must be defined for each calibration if you want to compare the measurements from one X-Ray pair to another camera or X-Ray pair. For example, the Event Summary document, that is described later, compares the results from multiple measurement documents to compute velocity and angular velocity. In order to compare these results, all the coordinates must be transformed into the same reference frame (the world reference frame).

Measurement

After a calibration has been completed and a test has been conducted. The resulting images from the test can be used to perform measurements.

A Measurement Document is created by selecting $File \rightarrow New$. Then select **Measurement**.

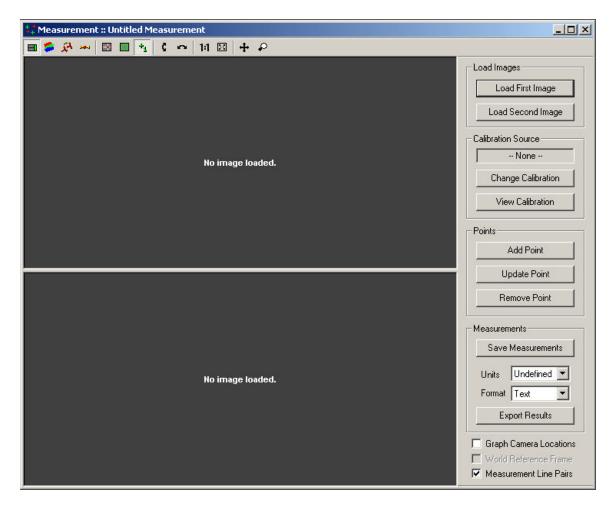


All the controls are located on the toolbar or on the panel on the right. The toolbar controls were described earlier in this document.

A pair of images from an actual test should be loaded into the Measurement document. An image from the top X-Ray system should be loaded using the **Load First Image** button. An image from the side X-Ray system should be loaded using the **Load Second Image** button.

To perform measurements, a calibration must have already been performed for the X-Ray systems being used. If a default calibration had been set, the default calibration is automatically loaded. If a default has not been set, then a calibration must be loaded by using the **Change Calibration** button.

The units of measure and the world reference frame transformation are also loaded from the calibration file. The units of measure that are displayed can be changed by changing the units in the panel on the right. The world reference frame transformation can be viewed or modified using the toolbar button or the right-click context menu.

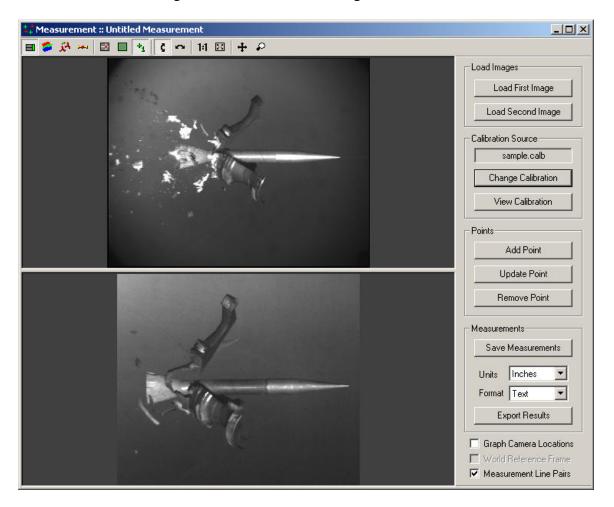


If you would like to review the calibration information, select the **View Calibration** button.

Measurements can either be the 3D location of a single point, or the 3D locations of the end points of a *Line Pair*. A line pair is a grouping of two sequential measurement points (points 1 and 2, 3 and 4, 5 and 6, ...).

Each Measurement document can display information as single point measurements or as line pairs. To change the display, click on **Measurement Line Pairs** checkbox in the right hand panel. When the Measurement document is set to show information as line pairs, additional orientation and length information is computed for each line pair and displayed in the Points listing. In addition, the first point in a line pair is shown in green and the second point in a line pair is shown in magenta. The connecting line is shown in cyan.

Typical steps to perform a measurement:

- 1. Click on a feature point in the top image. Zoom in using the mouse wheel to accurately select a feature point. The mouse wheel will zoom in/out on the currently selected point. It is always a good idea to zoom in on a feature point to ensure that you have selected the proper point.
- 2. Click on the corresponding feature point in the bottom image. Zoom in using the mouse wheel to accurately select a feature point. Verify that the uncertainty in the resulting image coordinates is acceptable. The uncertainty is shown on the bottom of each view. The red, blue, and yellow points and lines should be near one another in both images.
- 3. If you are satisfied with the feature points and uncertainty, click on Add Point.
- 4. If you would like to update the coordinates of a point, select the point in the Points list. Click on the correct feature point in the top and bottom image, then click **Update Point**.
- 5. To remove a point, select the point in the Points list and then click on **Remove Point**.

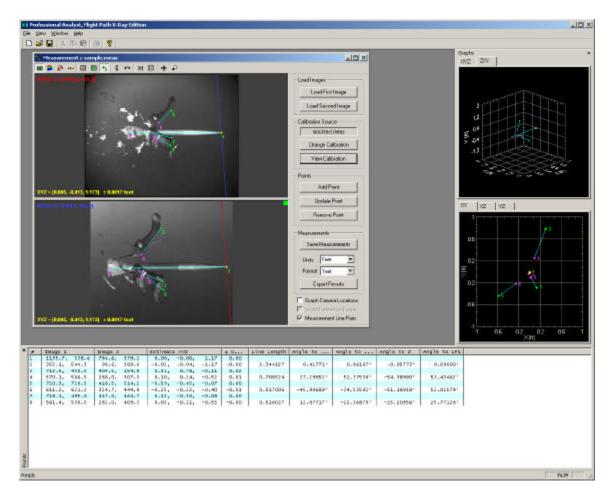

When you have selected all the points of interest, you can save your measurement results to a file. This will save the calibration information and measurement information to a measurement file (*.meas).

Note that the measurement file contains a copy of the calibration information. Once it is saved to the measurement file, the measurement does not use the original calibration file. If you make any changes to the original calibration file, you must reload the calibration into the measurement document using the **Change Calibration** button in order to apply the changes.

You can also export the measurement results to a Text document or to an Excel spreadsheet. First select the units and the export format and then hit the **Export Results** button. Here is a sample of the export results in Excel format. The sections are color coded to correspond to the first and second points in the line pairs and the line information for each pair.

3		yniat Iools				A1 21 40			line.		ten in 10				-	قلم .
A1.		Profession						· · U) .	Artal		10 10	B Z U		EB 1#	L • 💁 •	· 🔺 ·
A		C			F	G	H	1	L	к	L	M	N	0	P	Q
Profe	ssiona	Anal	vst. F	light												
-	Contentie		, out	S		. naj										
est Data		_	_	_	_	_	_						_	_		
ile Name:	D:\Samples	sample mea	20		_											
late:	5/28/2004	E														
	3:25:56 PM	-														
nits:	Feet															
leasuree	sent Data															
irst Point			_		_				Second Po		_			_	_	
nage 1		Image 2		Estimated	Pasition		± Uncertain	ntry.			Image 2		Estimated	Pasition		± Unce
			V .			Z		- C				V.			Z	10000
														-		1000
1135.65		784.556		0.001158					385.093 570.149	544.311 506.53			0.095806	-0.03942	-1.16821	
720.323				-0.58553					611.163					-0.22425		
718.148	495.848	416.953	464.734	0.132638	-0.29899	-0.081	0.001893		561.444	530.47	282.009	409.257	0.025906	-0.10686	-0.54788	-0.00
						_										
																_
		Sheet3 /														

If the same calculations are always done in Excel, the export feature can also be told to use an existing Excel file as a template for the export results. This template can contain formulas, graphs, or any other Excel elements. Select **Excel Template** as the **Format** type before exporting. Then select the Excel file to use as a template.

Event Summary

After you have created a set of Measurement documents for one event, you can combine these all in one Event Summary Document. An Event Summary document loads a set of Measurement documents and displays a summary of all the measurements contained in each of the Measurement documents.

An Event Summary Document is created by selecting File \rightarrow New. Then select Event Summary.

µ∕ Event	Summary ::	Untitled Summary					_ 🗆 ×
	Time (sec)	File Name					
Meas 1				Load	Edit	Reload	Remove
Meas 2				Load	Edit	Reload	Remove
Meas 3		_		Load	Edit	Reload	Remove
Meas 4				Load	Edit	Reload	Remove
Meas 5				Load	Edit	Reload	Remove
Meas 6				Load	Edit	Reload	Remove
Save	Display	Length 💌	Units Undefined 💌	Form	at Text	•	Export
J							

Load each measurement document by clicking on the corresponding **Load** button. For each measurement document, you must specify the time associated with that measurement. This allows the application to compute velocity information between measurements.

Up to six (6) measurement documents can be combined into an event summary. After you have loaded all the measurement documents, you can save the event summary information by clicking on the **Save** button. All the time and measurement document information will be saved to an event summary file (*.esum).

If you wish to make changes to a measurement document, you can hit the **Edit** button next to the measurement document file name. After you have finished making changes and have saved the measurement document, you must hit the **Reload** button so that the new changes are loaded by the event summary.

The Points listing will display information about each line pair contained in each measurement document. Which quantity is displayed is controlled by the **Display** drop down control in the Event Summary document window.

	98						_	Oats			
								Oats			
								XY2	28/ 1		
										TT	
-		NAME OF TAXABLE PARTY.	_					182.0	E-++		
	Event Summary 1							- 80.2		140	
	Time (sect	File Nane D VSaroleri vanple 1 wess	Load Eet	Reload R	arove						
	Hear? Ett	D Vserplezkoanple 2 teau	Load Idt	A resourced into	CHUNE .			178.6			
	Heat 3 E.O2	D Sampler/cample 3 wear	Load Edt		erove			-			24.1
	Meas 4 1.03	D VSamples/vample 3 wear	Load the		engye				A		100 C
	Head 5 ED4	D Viseples/sample 5 secs	Lowd Eek	a success of the	07010						
	A DOWNER WATCHING TO A DOWNE WATCHING TO A DOWNER WATCHING TO A DOWNE WATCHING TO A DOWNER WATCHING TO A DOWNER WATCHING TO A DOWNER WATCHING TO A DOWNER WATCHING TO A DOWNE WATCHING		and the second second	annexed an							
	Mean 6 LUS	D. \Semples\cample 6. seec	Lond Edit	Reloat R							
			a de la constante de la constan	1	enove			W	12 12	1	
	name i anno	(20V 718	-	1	
	Sine Diploy	Foster 🗶 Bole Inches	Fanal Fed		tapos				-		
	Save Diploy	Foster 🔳 Bolts Inches	Fanal Fed					716			
	Sine Diploy	Frailan 🕑 Ueits Inches	T Fanal Ted								
	Sine Diploy	Faster 🗶 Uols (Index	■ Fanal [red					223.6			
	Sana Dipite	Faster 🗶 Uols (Index	Fanal Fee					716			
	3 me Nobe	Foster 🔳 liek Index	▼ Fand [Fed					718 3238 6712 32			
	Sine Diploy	Foster 🔳 Unit Index	T Fanat Fed					718 3238 4712			
	Sine Diploy	Fradien 💌 Ukás Jindes	Tanki Fer	2				716 3336 471 2 3488 3488			
ŀ	Sire Digity	Frater 🕊 Uras Inches	Fanal Fee					718 3238 6712 32			
	3aw Diple	(Foster 🗶 Urde Jinches	Tons Ter					778 933.6 471.2 5 5 349.8 228.4		•	
	Save Digits	(Foster 🔳 Units (Inches	Tona Ter					778 2236 671 2 3 3 9 9 9 9 9 9 9 9 9 9 9 9 9 9 9 9 9			175.5
	Save Digits	Fradien 💌 Ukás (Inches	Tonal Fer					778 2236 671 2 3 3 9 9 9 9 9 9 9 9 9 9 9 9 9 9 9 9 9		•	175.6
6 March 1	70.4		76.05			964		716 93336 671 2 36 5 3488 235.4 104	BIL AIR6	61 2 57.2 2000 [a]	173 6
-2,74, 4,85	». 104.17 -5.	5 Prior 7 7 4.75 204.4 - 75.75 4.		4. 5, 434	Mass 1 4 -3.7, 4	.2, 504.4 -1.	4.7.	716 2236 6712 3.68 235.4 104 2704.4	nil Air6	61 2 57.2 200 [a]	173 6
-2.74, 4,85 -2.58, 6,42 -5.28, 4,89	7, 104.07 - 7, 1, 105.66 - 2, 1, 105.66 - 2, 105.10 - 100 - 2,	5 1 Priori 7 7, 4, 5, 204,45, 7, 4, 6, 6, 4, 205,92, 6, 6, 4, 10, 205,92, 6, 6,	96-46 5 204-4 - 5,7, 4 205-9 -2;6, 1 205-9 -2;6, -1 -4, -2 -4, -2 -4, -2 -4, -4, -2 -4, -4, -4, -4, -4, -4, -4, -4,	4. 5, 434 6. 4, 435 8. 0, 435	Maas 1 4 -3,7, 4 9 -3,-5, 6 9 -1,-5, 6	-5, 504,4 -1, -4, 506,9 -3, -0, 809,3 -5.	4,5 6,4 1.0,	776 2236 471 2 35 3466 275 4 104 2 704, 4 705, 9 705, 9	still Attros	St 2 Milli Second	
-2.58, 6.42 -3.28, 4.88 -1.78, 3.52	P. 104-1753. 5. 005.06 - 2- 7. 008.095. 105.87 - 5.	B PRIME P 7, 4, 55, 204, 4 75, 75, 44, 55, 75, 74, 74, 75, 75, 75, 75, 75, 75, 75, 75, 75, 75	Model 15 2034.4 -5.7 14 207.9 -5.7 15 2037.9 -41.6 15 207.9 -5.7	4.7, 434 6.4, 455 8.0, 485 3.7, 485	Mass 1 4 - 2-7,7 4 5 - 2-7,6 6 9 - 2-7,7 9 6 - 3,7,7 9	-5, 504,4 -1, -4, 506,9 -1, -0, 809,9 -5, -5, 508,8 -1,	4,3, 6,4, 8.0, 3,3,	776 2236 277 2 5 5 5 3488 2764,4 704,4 704,4 704,4 704,4 704,4 704,4 704,4 704,4 704,4 704,4 704,4 704,4 704,4 705,5 705,8	HI ARG	61 2 57.2 200 [k] indies indies	
-2,74, 4,85 -2,58, 6,42 -6,28, 4,99 -1,78, 3,52 -1,42, 5,82 -1,68, 2,23	7, 104.07 - 7, 1, 105.66 - 2, 1, 105.66 - 2, 105.10 - 100 - 2,	2 PROF 2 7 4.5 204.6 3.7, 4. 4. 5. 204.6 3.7, 4. 4. 5. 201.9 2.7, 5. 5. 5. 5. 501.8 5.7, 5. 5. 5. 6. 201.8 5.7, 5.	96-46 5 204-4 - 5,7, 4 205-9 -2;6, 1 205-9 -2;6, -1 -4, -2 -4, -2 -4, -2 -4, -4, -2 -4, -4, -4, -4, -4, -4, -4, -4,	4.5, 434 6.4, 435 8.0, 440 5.5, 434 5.5, 434 2.1, 440	Nage 1 4 - 2,7,7 4 4 - 2,7,7 4 5 - 2,6,7 5 7 - 2,7,7 5 9 - 2,7,7 5	-5, 504,4 -1, -4, 506,9 -3, -0, 809,3 -5.	4,5, 6,4, 8,0, 3,3, 5,0, 2,2,	7716 223.6 671 2 5 7 3.468 23.6 3.468 23.6 4 7 0 4.6 7 704.4 7 705.8 704.5	hit Altr6	66 2 57 2 200 [bi] Units Units Units Units	

The Event Summary results can also be exported to a Text document or to an Excel Spreadsheet.

Select the units and the export format and then hit the **Export** button. This will export all available information about the measurement line pairs. The current display setting is not used during export. All possible display values are exported simultaneously.

	CALL COLORS	icel-Bookt Vew (rset Fo	and Tasks	Data Useda	AL LOUIS A												کلم فلم
-	a the state of the local data		and a local plant of the local plant		the second s		AL 21 4	tars	. 🗊	and .		10				. d.	
				nal Analyst					7.000	1 com		110 218-	4 2		100 No.	H	-
	A	8	Professio	D D	E	F	G	H	1		к	1		N	0	P	Q
i												-	18	19	0		
	Profe	ssiona	al Ana	lyst, F	light	Path	X-Ray	Editio	on Da	ta She	et						
	0 B	1-2 X	1	100 200	S.S.C. 1	1			-	1	-	1		-	1	1.	
	con sub																
1	Test Data																
		12410 200															
		DOSamples		um													
	Date:	5/28/2004															
	Time:	3:24:03 PM															-
	Units:	Inches															
	deasures	nent Files															
	Vieas 1	Time	0	File Name:	0.\Samele	sisample 1	meas										
	Meas 2	Time		File Name:													
	Meas 3	Time:		File Name:													
1	Meas 4	Time	0.03	File Name:	D:\Sample	sistemple 4	meas										
	Weas 5	Time		File Name:													
l	Meas 6	Time	0.05	File Name:	D:\Sample	is/sample 6	meas										
		A PLO															
Ļ	10000																-
	ength	Sec. 1	Marca 7	Marso 7	Maren K	Marca F.	Marco D										-
1	ine	Meas 1 11.523129			Meas 4	Meas 5	Meas 5										
ř				5.002484													-
		5.0037875															
		4.9964411															
		3.1109962															
	E	2.4222407	2.422241	2.422241	2.422241	2.422241	2.422241										
	7	3.2668137	3.266814	3.266814	3.266814	3.266814	3.266814										
	Position																
L	.ine			Meas 1 Z													
		-3.7367562															
		2 -2.5826431				6.421658		-2.58264				6.421658	405.856		6.421658		
		-3.6971304				3.319653			3.319653			4.907000			4.907000		-0.38
		-5.4156496															
		-3.6890189						-3.68502								504.8634	
		-2.8023769														504 8802	
l	Displacen	nent						1									
	.ine	Meas 1			Meas 4	Meas 5	Meas 6										
		104.5486				504.4847											
		105.08209															
		106.15811															
		105.95072															
		105.17136		305.0011	404.9791	504.965/	704,9604		-	·							-
	F N.Sh	est1 / Sheet2	/ Sheet3 /								4					Contraction of the	Concession in which the

If the same calculations are always done in Excel, the export feature can also be told to use an existing Excel file as a template for the export results. This template can contain formulas, graphs, or any other Excel elements. Select **Excel Template** as the **Format** type before exporting. Then select the Excel file to use as a template.

	icel - Bookt															als.
and the state of the local data	Vew Insert R	and a local production of the local of	and the second second second	the second se			-									فلم
	3 8 4 3							· · [2] •	Artal		16 - 1	a Y U		田保	田• 3.	· ≙ ·
A1			nal Analyst													
A	В	C	D	E	F	G	н	- Arrest		К	L	M	N	0	Р	0
Profe	ssiona	al Ana	lyst, F	light	Path 2	X-Rav	Editio	on Dat	ta She	et						
-	TA S			1000			C				_			_		_
Test Data	1															
10.00	2012	100														
File Name	a:D:\Samples		um													
Date:	6/1/2004															
Time:	4:56:00 PM															
Units:	Inches															
	and Piles															
Measuren Meas 1	nent Files Time:	0	File Name	0.)Samel	ala anala 4	Print Part										
Meas 2	Time		File Name						-	_				-		
Meas 3	Time		File Name								Line1				-	-
Meas 4	Time		File Name												-	
Meas 5	Time		File Name:					800 -								
Meas 6	Time		File Name					700			_		-			
	1.000			-14100-1417	100.00			600	0			1.10			-	
								8 900 400	-			1	-		-	
Length								400			//	<u></u>	28	Position		
Line	Meas 1			Meas 4	Meas 5	Meas 5		300			/	-		placement		
	11.523129							E 200		//					-	
3			5.002484					100	1	~					-	
			5.003788					100	/						-	
	4 4.9964411 5 3.1109982		3.110998					32	1 2		4 8	6 7	2.0		-	
	5 2.4222407								- S		roment	CO 25				
	3.2668137									Principal D	Contraction of the local distance of the loc				-	
	0.2000104	0.200014	0.200014	3.200014	3.200014	3.200014		<u> </u>								
Position																
Line	Meas 1 X	Meas 1 Y	Meas 1 Z	Meas 2 X	Meas 2 Y	Meas 2 Z	Meas 3 X	Meas 3 Y	Meas 3 Z	Meas 4 X	Meas 4 Y	Meas 4 Z	Meas 5 X	Meas 5 Y	Meas 5 Z	Meas B
	-3.7367562															
	2 -2.5826431	6.421658	105.856	-2.58264	6.421658	205.856	-2.58264	6.421658	305.856	-2.58264	6.421658	405.856	-2.58264	6.421658	505.856	-2.58
	3 -5.3910453										4.987868				505.9017	
	4 -3.6971304				3.319653			3.319653			3.319553			3.319553		-3.69
	5 -5.4156496														504.9117	
	3.6850189					204.8634			304.8634		3.327607				504.8634	-3.68
7	-2.8023769	6.405605	104,8602	-2.60238	6.405605	204,8802	-2.60238	6.405805	304 3602	-2.60238	6.409805	404,8602	-2.60238	6.409805	504,8802	-2.80
Disalar	Meas 1	Meas 2	Meas 3	Meas 4	Meas 5	Meas 6										
			304.4294		504.4847											
	104.5498															
Line 1		205 9724				100,00										
	2 105.08209					706 9999										
Line 1	2 105.08209 3 106.15611	206.0327	305.9344 305.9899 305.8745	405.9682	606.9651											
Line 1 2 2	2 105.08209 3 106.15611	206.0327 205.8941	305.9899 305.8745	405.9682	505.9651 505.8585	705.8516										

All the information is exported into the template file in Excel, with any calculations or graphs being automatically populated.